



Software installation guide

Step 1, ECS application installation


Open CD, Run ECS10SETTUP.EXE:

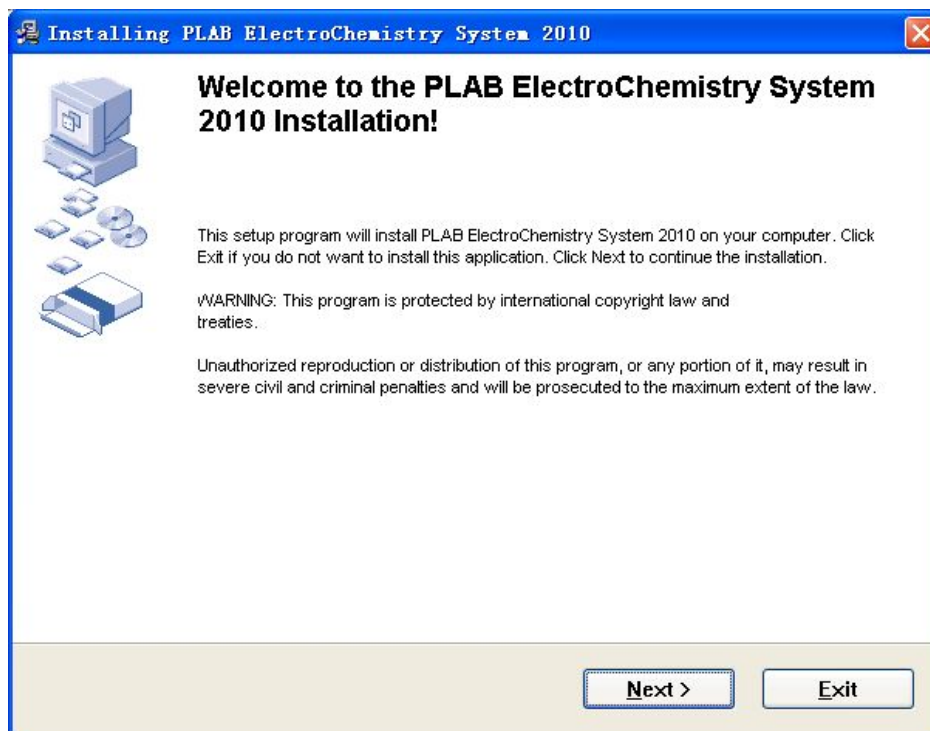


Click 'Next'

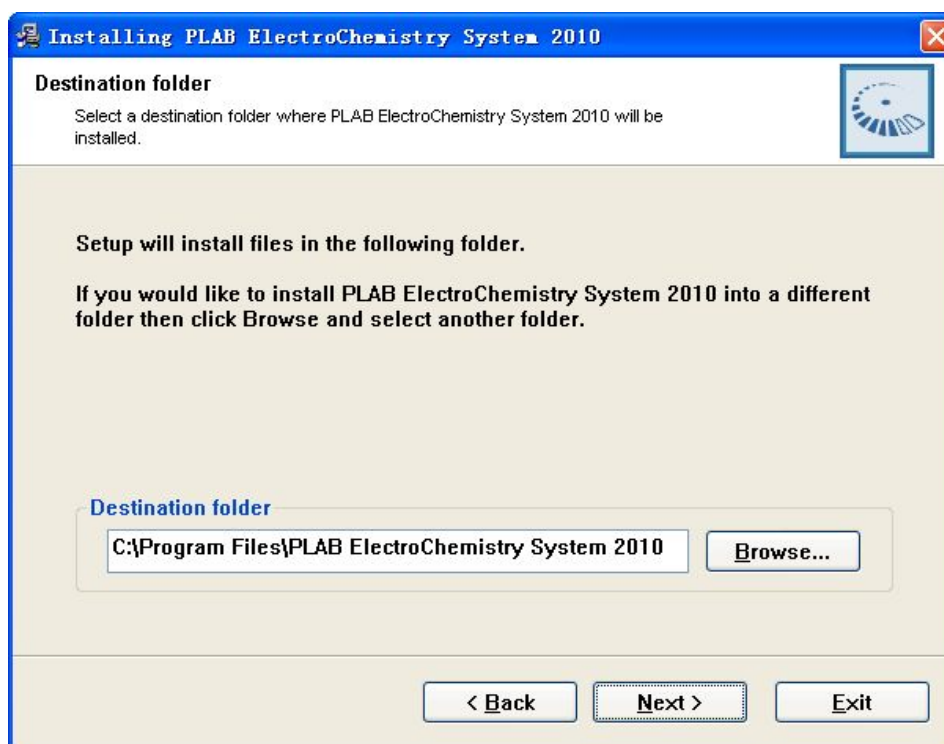

If the following window appears, click 'Overwrite All'

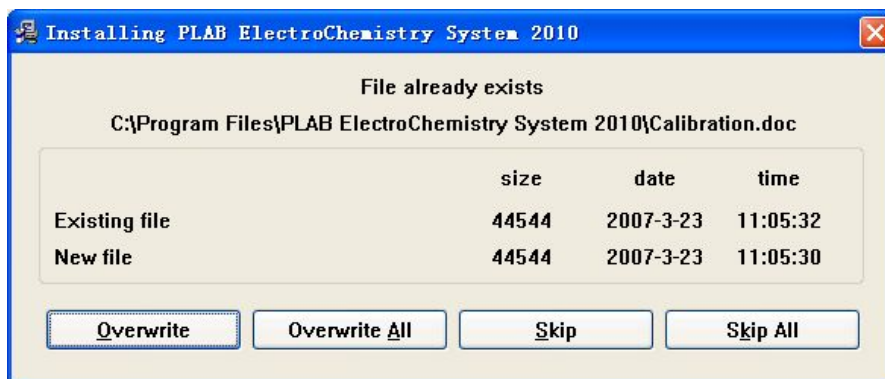


Click 'Finish' to finish the application installation and driver installation will start automatically.

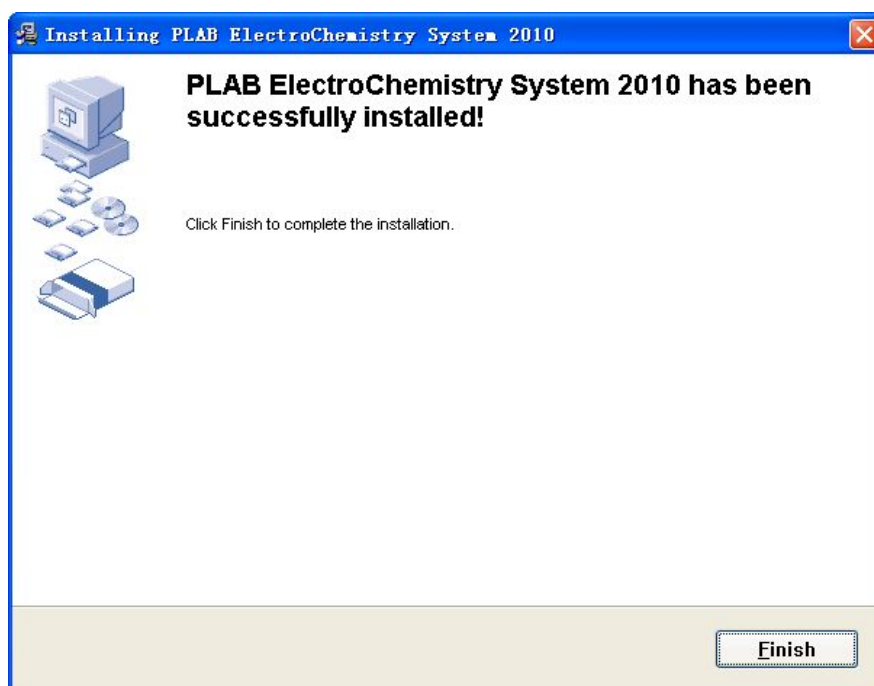


Step 2, Base driver installation

Switch to driver installation window, press any key to start

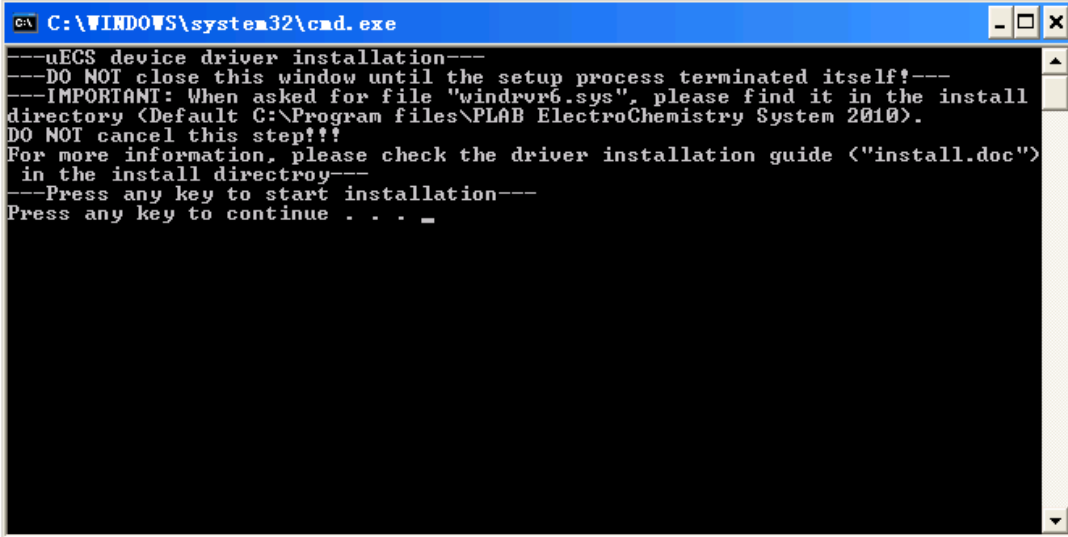

```
C:\WINDOWS\system32\cmd.exe
--uECS device driver installation--
--DO NOT close this window until the setup process terminated itself!--
--IMPORTANT: When asked for file "windrvr6.sys", please find it in the install
directory (Default C:\Program files\PLAB ElectroChemistry System 2010).
DO NOT cancel this step!!
For more information, please check the driver installation guide ("install.doc")
in the install directroy--
--Press any key to start installation--
Press any key to continue . . . _
```

If the following prompt appears, unplug uECS , then press 'R':

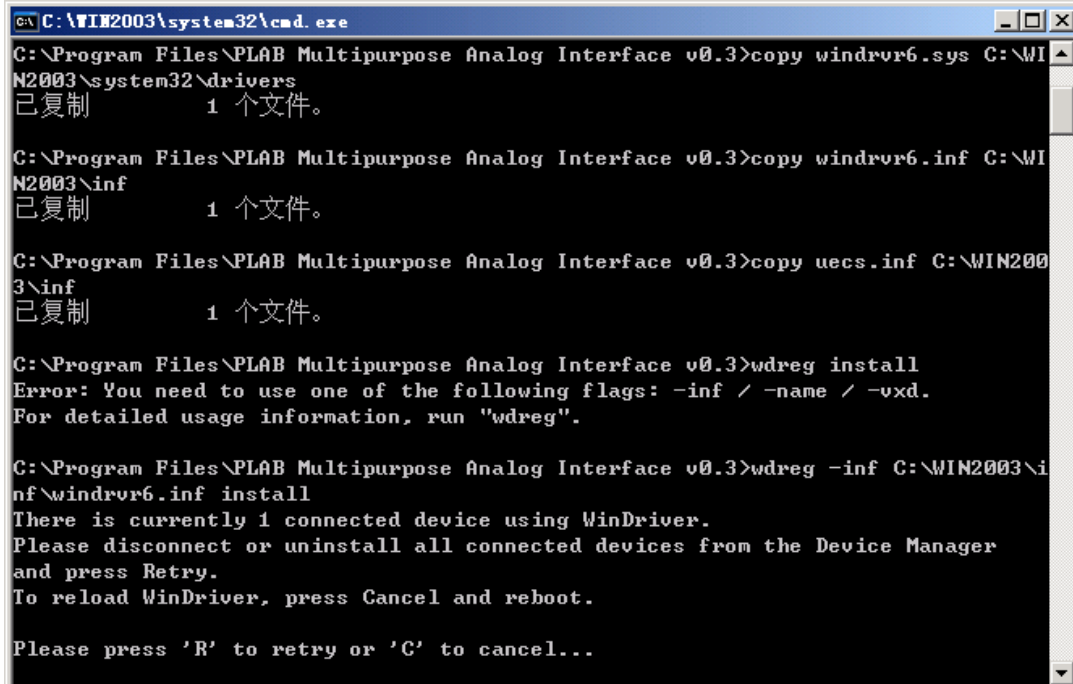

```
C:\WIN2003\system32\cmd.exe
C:\Program Files\PLAB Multipurpose Analog Interface v0.3>copy windrvr6.sys C:\WI
N2003\system32\drivers
已复制      1 个文件。

C:\Program Files\PLAB Multipurpose Analog Interface v0.3>copy windrvr6.inf C:\WI
N2003\inf
已复制      1 个文件。


C:\Program Files\PLAB Multipurpose Analog Interface v0.3>copy uecs.inf C:\WIN200
3\inf
已复制      1 个文件。

C:\Program Files\PLAB Multipurpose Analog Interface v0.3>wdreg install
Error: You need to use one of the following flags: -inf / -name / -vxd.
For detailed usage information, run "wdreg".


C:\Program Files\PLAB Multipurpose Analog Interface v0.3>wdreg -inf C:\WIN2003\i
nf\windrvr6.inf install
There is currently 1 connected device using WinDriver.
Please disconnect or uninstall all connected devices from the Device Manager
and press Retry.
To reload WinDriver, press Cancel and reboot.

Please press 'R' to retry or 'C' to cancel...
```


If the following window appears, press “Yes” (Red circled button)

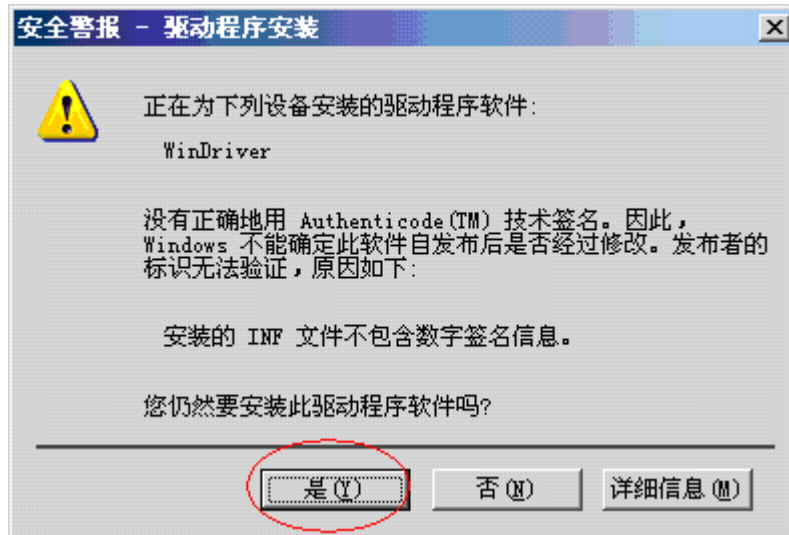


When this window appears, press "Yes" (Red circled button)

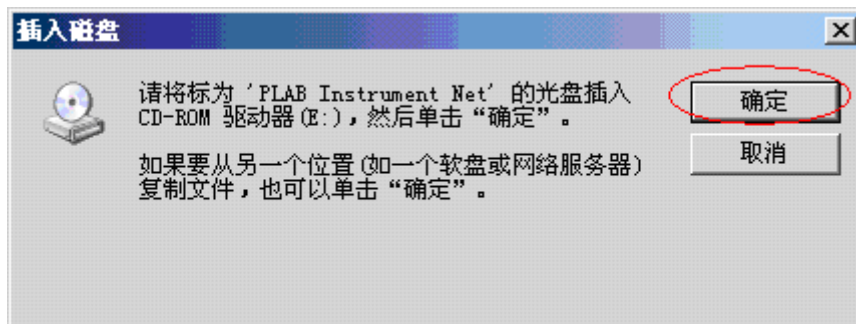


Press "Browse" (Red circled button)

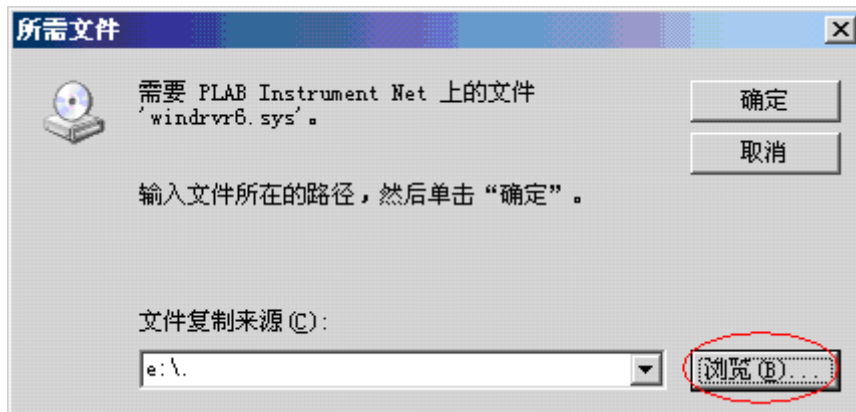

Select the installation directory (Default C:\Program Files\PLAB ElectroChemistry System 2010), then press "Open" (Red circled button)

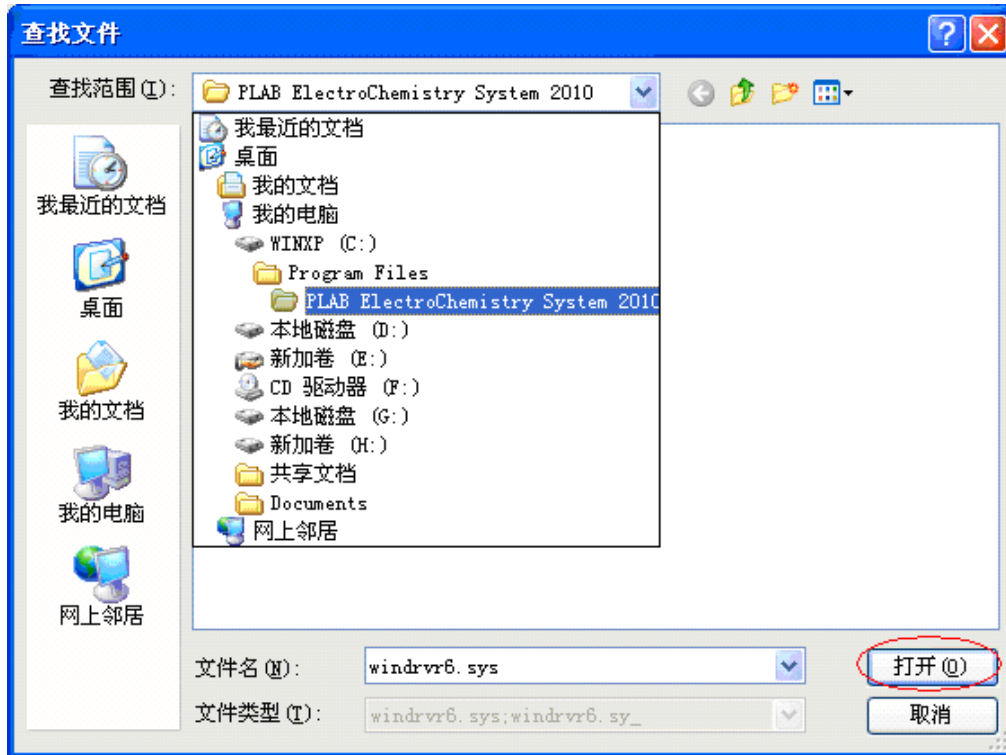


Press "Yes" (Red circled button), now base level driver is installed.

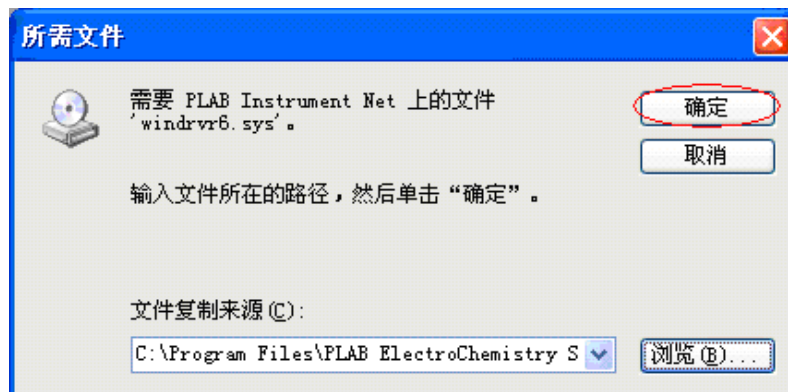

Step 3, uECS driver installation


When uECS is plug in a USB port for the first time, the following window will appear, select “Automatic installation driver”(the first option), then press “Next” (Red circled button)

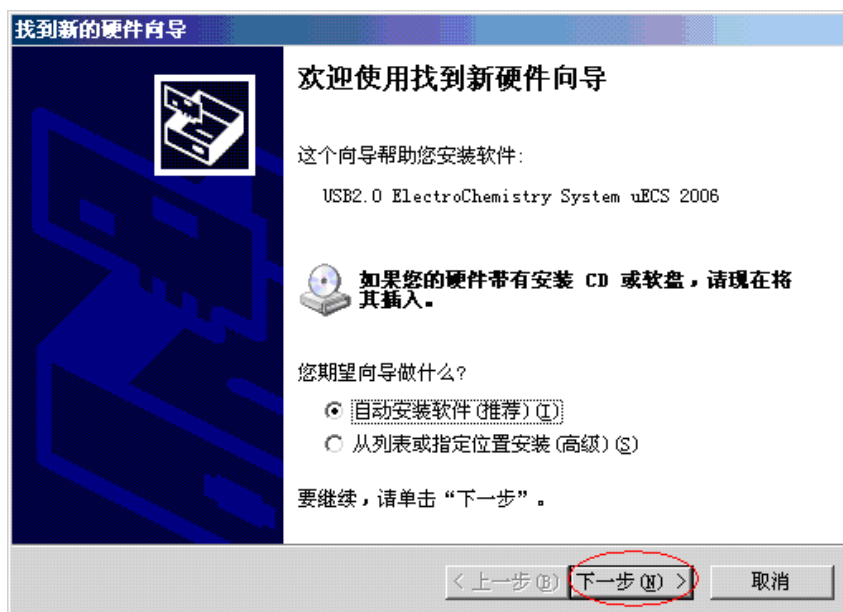

Wait



Press “Yes” (Red circled button)

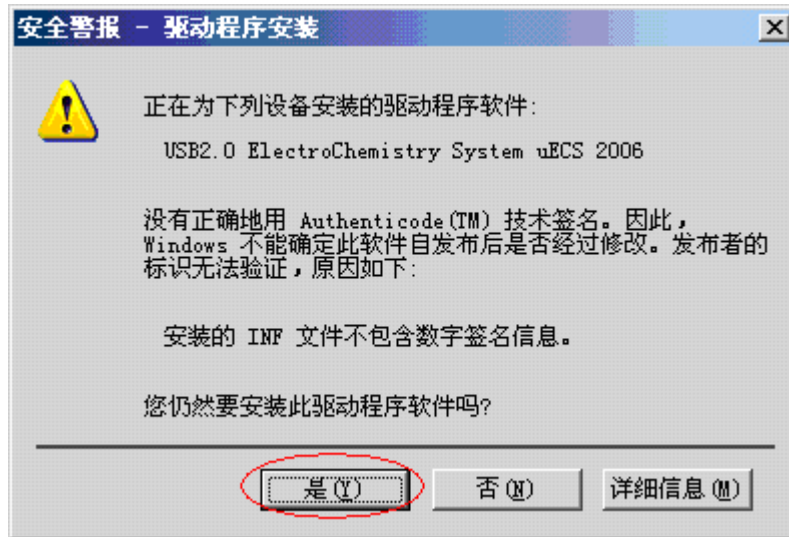

Press "Finish" (Red circled button)



Step 4, Checking driver installation

Right click "My computer", select "Manage"(high-lighted menu item).



Click "Device manager" in left side.




Browse "Jungo" folder by pressing "+" button



If both “USB2.0 ElectroChemistry System uECS 2010” and “WinDriver” appears without exclamation marks, it means the driver installed successfully.


Step 5, Checking and updating License

Start ECS program, a splash window including license information will appear for a few seconds. It can also be open by clicking “Help-About” menu item:

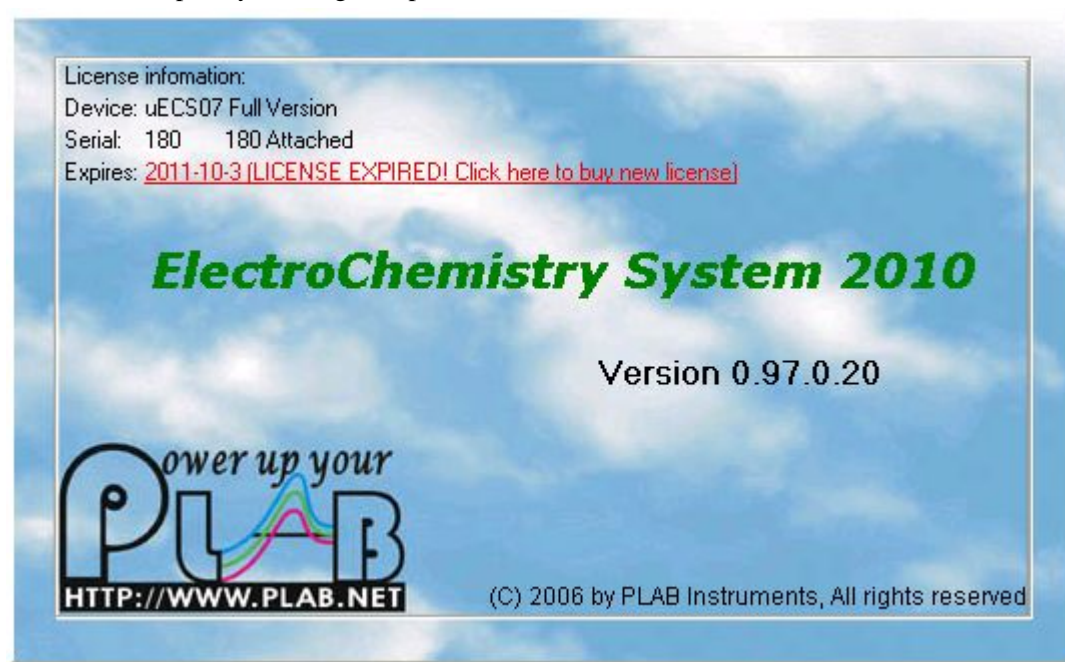


The first line of license information is the device type of current license code. The second line is the serial number of your current license code and the serial number of the attached device. The 3rd line shows the date when your license expires.

If the current license code is expired, you should get a new code from the provider, then update license code by pressing “Help-UpdateLicense” menu item:

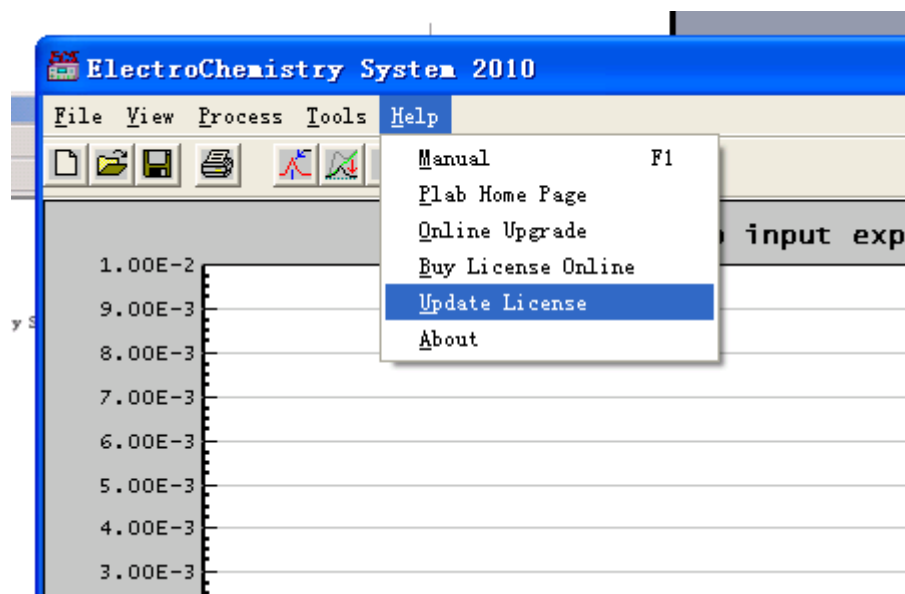

Input new license code in this dialogue and press OK:

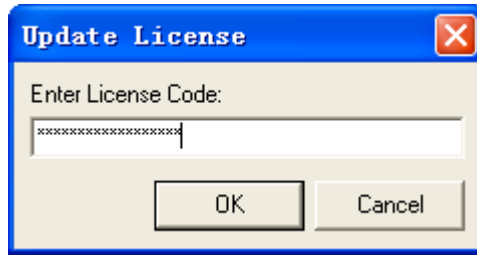

After updating license code, restart ECS program and check the license information in the splash window.